

Poly-L-Lysine Coating of Coverslips

1. Dilute Poly-L-Lysine solution 1:10. Aliquot 300 μ Ls of poly-l-lysine solution onto coverslips (enough to cover surface of coverslip).
2. Incubate coverslips in poly-l-lysine solution for 25 minutes, under sterile hood.
3. Aspirate solution from coverslips, and allow to air-dry for 25 minutes under sterile hood.
4. Use coverslips immediately, or cover and store in dark, dry space for 1 week before use.

Materials:

- Poly-L-Lysine Solution (Sigma Aldrich, cat # P8920)